






GW4260 – GW4290 INSTRUCTION SHEET

SELECTING AND STARTING A WASHING PROGRAM



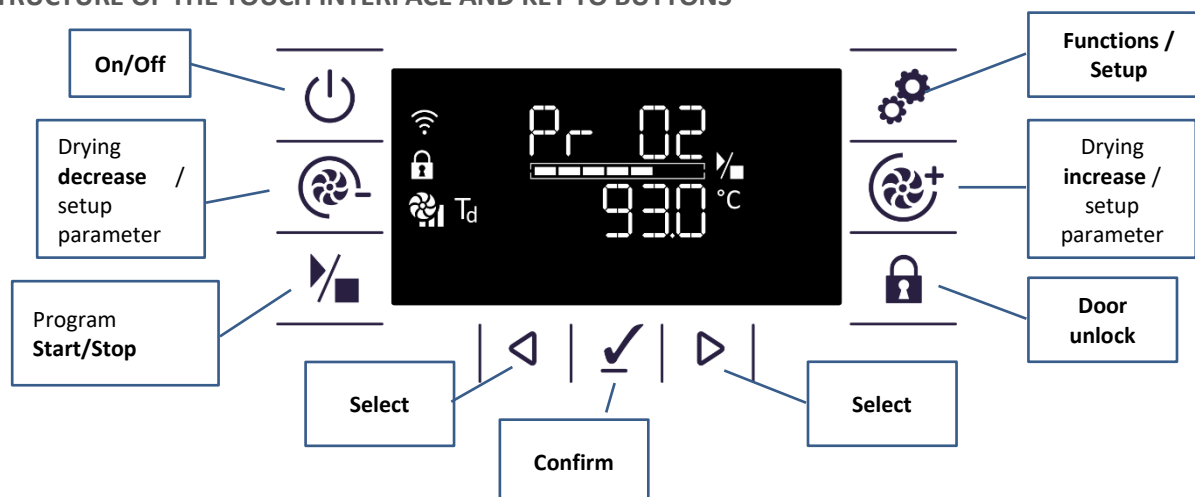

	Switch on the device by pressing the On/Off button (hold down for 2 seconds). If the door is open: close the door by shutting it against the device until the lock clicks shut. The buttons are only enabled with the door closed. [For GW4290 only: power up the device by turning the main switch from “0” to “1”. The switch is in the side cabinet.]
	Select the required program with the Selection buttons.
	Press the Start/Stop button – hold down (2 seconds).
	Only once the program has been successfully completed: the second line of the display shows “End”. Unlock the door with the “padlock” button to open the door, extract the trolley and collect the load.

STRUCTURE OF THE TOUCH INTERFACE AND KEY TO BUTTONS



N.B.: in models without drying system, the associated buttons increase and decrease the setup parameters; the icons are differentiated.

SPECIFIC ALARMS AND WARNINGS

ICON	MEANING
	Alarms are signalled by means of the icon with the “Danger” symbol and a code on the display. “AF” – fatal alarm. “W” – Warning, non-fatal alarm. Press Confirm to acknowledge an alarm or warning. When an alarm occurs, the device responds to the problem itself. The number of the warning or alarm appears beside the letter code: refer to the manual for correct management of the event.
	Low detergent level: replace the empty jerry can with a new one.
	Machine maintenance: routine maintenance required.
	Service Icon: indicates that program is for service and not for load processing.

PROGRAMS

1 Prewash	11 GW Thermal disinfection 93 °C – 10' + Drying
2 Plastics 75 °C – 3' + Drying	12 Intensive GW thermal disinfection 93 °C – 5' + Drying
3 Fast Glassware 75 °C – 3' + Drying	13 Intensive GW thermal disinfection 93 °C – 10' + Drying
4 Medium Glassware 75 °C – 5' + Drying	14 GW Petroleum Dirt 93 °C – 1' + Drying
5 Standard Glassware 80 °C – 5' + Drying	15 Intensive GW Petroleum Dirt 93 °C – 5' + Drying
6 Intensive Glassware 90 °C – 3' + Drying	16 GW Petroleum Grease 93 °C – 1' + Drying
7 High temperature glassware 93 °C – 3' + Drying	17 GW Mineral Oil 93 °C – 10' + Drying
8 Glassware Agar 93 °C – 3' + Drying	18 GW Petrol 93 °C – 5' + Drying
9 Intensive Agar 93 °C – 5' + Drying	19 GW Diesel 93 °C – 1' + Drying
10 GW thermal disinfection 93 °C – 3' + Drying	20 GW Universal Petrol Diesel 93 °C – 10' + Drying

THIS DOCUMENT

Brief instruction sheet for basic procedures. For details relating to correct, safe use of devices, the updated program data, interpretation of alarm messages and installation instructions: refer to the full version of the user manual. **N.B.:** the device must be commissioned by authorised technical staff; the recommended detergents and dishwasher salt are required for this purpose. For periods out of use: consult the manual to prevent problems in the device such as crystallisation of liquid detergents.